PREVENTATIVE MAINTENANCE CHECKLIST FOR DIRECT HEAT AND WATER JACKET INCUBATORS

MfgModel/Serial	The server of the least
Customer	Thermo Fisher
Location	
Contact / Phone	•
Contract Billable Oth	
Complete all applicable tasks with the following codes and comments. Include additional comments on rear. P Pass F Fail N N/A	
TASK	COMMENTS
Inspect	
Inner Door Gasket	
Outer Door Gasket	
Inner Door Latch	_
Remove / Clean	
Humidity Water Pan	
Interior Shelves	
Chamber HEPA	
Ducts and Plenums	
Air Sample Filter Blower Motor Assembly and Gaske	<u> </u>
CO2 Sensor and Gasket	t
CO2 Filter	
Access Port Filter	
Inner Door Gasket	
Using a non-chlorine based genera	l laboratory cleaner, wipe out entire chamber, including door
opening, inner door gasket, plenum	, ducts, shelves, blower cavity and sensor cavity
Rinse twice with clean water	
Dry the entire chamber	
Landall .	
Install CO2 Sensor and New Cocket	
CO2 Sensor and New Gasket New Blower Wheel	
Blower Assembly and New Gasket	
New Access Port Filter	
New CO2 Filter	
Inner Door Gasket	
Plenum and Ducts	
New Air Sample Filter	
New Chamber HEPA Filter	
Shelves	
Water Pan with Fresh Water	
Inner Door Gasket	
Rust Inhibitor (WJ only - every 2 ye	ars) nimum of 8 hours and verify all parameters.
	minum of 6 hours and verify all parameters.
-	
Verify	
Over Temperature Alarm	
Low Temperature Alarm	
High CO2 Alarm	
Low CO2 Alarm	
Door Ajar Alarm Temperature Sensor Failure Alarm	
CO2 Sensor Failure Alarm	
Low Water Alarm (WJ only)	
Type of Water / Additives	
Engineer:	Date: